

(continued) Appendix B. Repairer PMCS

4110-01-287-7111
Refrigerator, Solid State, Biological, Model RCB42P

[M-Monthly, Q-Quarterly, S-Semiannually, and A-Annually]

ITEM NO	INTERVAL	ITEM TO BE INSPECTED AND PROCEDURE	IS NOT MISSION CAPABLE IF:
1	S	<p>Refrigerator, Solid State, Biological</p> <p>Conduct an inventory to ensure that the items listed on the Equipment Parts and Accessories List are on hand.</p>	<p>Missing components or accessories prevent the operation of the refrigerator.</p>
2	S	<p>Maintaining Refrigerator</p> <p>a. Verify that the refrigerator functions as directed by the Operators Preventative Maintenance Checks and Services manual.</p> <p>b. Verify that the unit functions on AC current.</p> <p>c. Verify that the unit functions on DC current.</p> <p>d. Check screw connections as directed by the manufacturer's service manual.</p> <p>e. If necessary, conduct the "ACCU" as directed by the manufacturer's service manual.</p> <p>f. Verify the electrical safety.</p>	<p>The refrigerator cannot function on AC.</p> <p>The refrigerator cannot function on DC.</p> <p>The refrigerator fails any of the electrical safety tests.</p>